


Powerful, high-performance measurement solutions

Essential Potentiostats



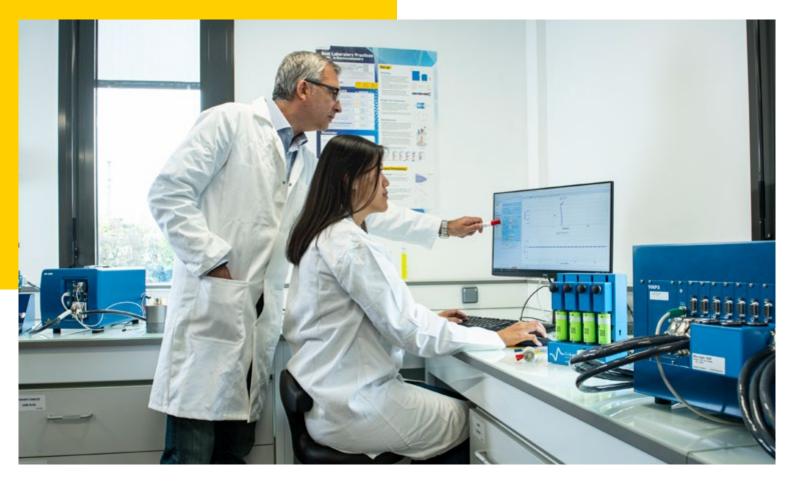

Benchmark software. High-performance hardware.

The perfect marriage of performance and control.

Modular, durable, single and multi-channel electrochemical workstations designed to meet both classical and demanding electrochemical research needs, the Essential range features EIS up to 1 MHz, a current range from 1 A down to 10 μ A and the possibility of extending to 800 A and down to 1 nA with modules.

No compromise has been made on quality in a range of potentiostats driven by EC-Lab® software, the same control interface that drives BioLogic's Premium instruments.

Advanced functionality such as Ethernet capability will help you manage multiple instruments from one computer as well as facilitate group-working. And **Quality Indicators** make the validation of EIS measurements simple.

Visit our YouTube channel and Learning Centre for scientific articles, EC-Lab® tutorials and product support information. https://www.biologic.net/topics/



Essential measurement tools

for electrochemists

Specification

Minimum/Maximum Capabilities

Channels: 1 to 16

Standard Voltage: ±10 V

Max Voltage: +60 V

Current: 1 nA to 800 A

EIS: Up to 1 MHz

Multichannel




Single channel



Ethernet capability for increased flexibility/improved group working

Share an instrument's channels via multiple PCs, or share multiple instruments from your PC.

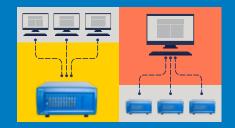

The most comprehensive and user-friendly software available

EC-Lab® is widely regarded as the **benchmark control and analysis software** by scientists across the globe. As simple to use as it is powerful, EC-Lab® offers a wide range of unique functionality that can help drive your projects forward.

Preset or bespoke techniques

80+ techniques

Integrated graphics

Customize graphs within EC-Lab®

Experiment sequence builder

Build sequential experiments based on conditional limits

External device control

(multiple devices)

Extensive analysis functionality

Including Z Fit for advanced EIS modeling

No need to plan experiments

Modify measurements "on the fly"

SP-50e: The perfect instrument

for education & general use

Channel capability: 1

Voltage: ±10 V

Current: ± 1 A down to 20 nA



- Highly cost-effective workstation- now with EIS
- Simple, easy-to-use, compact chassis
- Ideal for general use and education perfect for new researchers
- Exploit the power and performance of EC-Lab® with a limited budget

SP-150e: From μA to 800 A The power to do more

- Two channels: perfect for rotating ring disk electrodes (RRDE)
- Highly modular potentiostat. Users can easily add low/high current boosters (no factory upgrade required)
- Future proof: instrument easily upgraded with EIS and a new channel



Channel capability: 2

Voltage: ±10 V

Adjustable between -20 V to +20 V

Current: ± 1 A down to 20 nA

VSP: A versatile, 5-channel, research-grade, instrument



- Versatile, modular instrument for general needs
- · Future-proof: instrument can be easily upgraded with EIS, low/high current options and new channels
- Optional 4 A booster available

VSP-3e: Tailor-made for energy applications

Channel capability: 8

Voltage: ±10 V

Adjustable between -20 V to +20 V

Current: ± 1 A down to 20 nA

 Compact, upright design reduces instrument footprint and saves valuable laboratory space

 Future-proof: instrument easily upgraded with EIS, low/high current options and new channels



VMP-3e: Versatility, power and performance. A do-it-all measurement tool.

Channel capability: 16

Voltage: ±10 V

Adjustable between -20 V to +20 V

Current: ± 1 A down to 20 nA



- Research-grade instrument with 16 channel capability
- Easily upgraded by user: add low-current modules
- Connect each potentiostat to an external high current booster channel perfect for battery research/testing
- Ethernet capability via LAN connect several computers/users to the same unit to facilitate group working


Add-ons: Customize your potentiostat to match your field of interest.

Modules by potentiostat

| Options | Specification | Application | SP-50e | SP-150e | VSP | VSP-3e | VMP-3e |
|-----------------------|---|--|------------|------------|------------|------------|------------|
| Low current option | LC option | Provides pA accuracy, for analytical electrochemistry, corrosion and also small battery cells | \otimes | \bigcirc | \bigcirc | \bigcirc | \bigcirc |
| High Power booster | 20 V boosters: 2, 5, 10, 20 A ±3 V at ±80 A 5 V at ±100 A 60 V at 50 A* 12 V at 200 A* *Up to 4 boosters can be connected to increase current capability | A Battery, supercapacitor, fuel cell, electroplating & electrolysis, Supercapacitor or fuel cell characterization Battery testing Battery pack characterization Large battery cells, supercapacitors or fuel cell characterization | \otimes | \bigcirc | ⊘ | \oslash | \Diamond |
| EIS | Up to 1 MHz | EIS measurements can be made and validated with BioLogic EIS Quality Indicators | \bigcirc | \bigcirc | \bigcirc | \bigcirc | \bigcirc |


Contour plots.

EIS

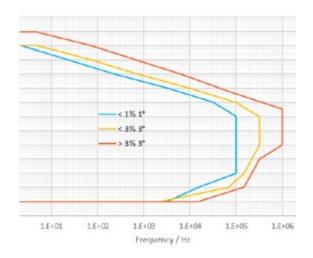

VMP3 contour plot (example applicable to rest of Essential Potentiostat range)

High Power

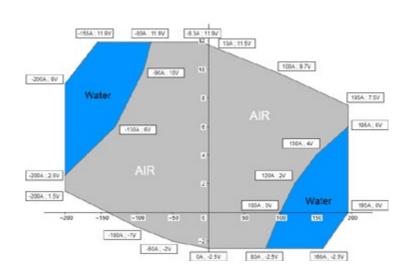

Contour plot demonstrating the high-power operating area of the FlexP boosters

Future-proofed potentiostats
Upgrade your own instrument, quickly, in your lab.
So your potentiostat grows with your needs.

Channel Specifications

| Features | | Essential Specification | | | |
|-----------------------|-------------------------|---|--|--|--|
| EIS capability | | 10 μHz to 1 MHz | | | |
| Analog Ramp Generator | | N/A | | | |
| Floating option | | CE to Ground | | | |
| Filters | | Software | | | |
| Acquisitio | n time | 20 μs | | | |
| Electrode | connections | 2, 3, 4, 5 | | | |
| IR comper | nsation | Manual, EIS, current interrupt (software) | | | |
| Current | | | | | |
| Maximum current | | ± 1A for "e-type" boards | | | |
| Current | With standard board | 6: 10 μA to 1 A | | | |
| ranges | With low current option | 9: 1 n A to 100 mA | | | |
| Lowest | With standard board | ±20 nA on 10 μA range | | | |
| accuracy | With low current option | ±2 pA on 1 nA range | | | |
| Lowest | With standard board | 0.8 nA on 10 μA range | | | |
| resolution | With low current option | 80 fA on 1 nA range | | | |
| Current | Internal | 4 A for VSP only | | | |
| booster | External | 2, 5, 10, 20, 80, 100 A, | | | |
| | | 200 A (FlexP0012), 50A (FlexP0060) | | | |
| Input impedance | | $1 T\Omega (//20 pF)$ | | | |
| Voltage | | | | | |
| Compliance | | ±10 V | | | |
| Max applied potential | | 0-20 V adjustable | | | |
| Resolution | | 5 μV on 200 mV | | | |
| Accuracy | | < 5 mV on ±2.5 V | | | |
| Maximum scan rate | | 200 V/s | | | |

Only with EC-Lab®

Modify-on-the-fly

No need to plan experiments
– you modify as you go, giving
you increased flexibility, easier
management of long-term
experiments and easier set-ups.

Full Cell Control

Measure (not only control) the voltage between positive and negative electrodes for batteries and fuel cells, just as you do with current.

Temperature Control Server

Manage climatic chambers from EC-Lab®, allowing users to perform automatic cycling with complex temperature profiles.

Z inst

Compensate drift during EIS measurements, for example, battery or specimens for corrosion studies.

Booster Specifications

| | 2///5.4 | 0/10/20 4 | 00 4 /UCD 007 | 100 4 /1100 1005 | FI - D00/0 | FL 00012 |
|---------------------------------------|---|--|---|--|--|--|
| | 2/4/5 A | 8/10/20 A | 80 A/HCP-803 | 100 A/HCP-1005 | FlexP0060 | FlexP0012 |
| Current | | | | | | |
| Compliance | 2 A: ±2 A, 4 A: ±4 A, 5 A: ±5 A | 10 A: ±10 A, 20 A: ±20 A | ±80 A | ±100 A | -50A; +49A | -200A ; +195A |
| Accuracy | | 10 A range: < 20 mA 20 A range: < 40 mA | | 100 A range: < 200 mA | 0.1%+/-0.01% FSR | 0.2%+/-0.02% FSR |
| Voltage | | | | | | |
| Compliance | Adjustable ±10 V range | Adjustable ±10 V range | ±3 V | 0.6 - 5 V | 0; 60V | -2.5; 11.9V |
| Control | ±20 V | ±20 V | ±3/5 V | ±3/5 V | 0; 60V | -2.5; 11.9V |
| Features | | | | | | |
| EIS frequencies | | 10 A: up to 80 kHz, 20 A: up to 80 kHz | up to 15 kHz, | up to 10 kHz | 10 kHz | 10 kHz |
| Floating mode | No | No | No | No | 5.6 kOhm | 5.6 kOhm |
| Rise/fall time (potentio, no load) | 15 μs | 25 to 60 μs | 95 μs | 1.7 ms | <10 µs | <20 µs |
| Parallel ability | No | _ No | No | No | Yes (up to 4 units) | Yes (up to 4 units) |
| Connection (terminal leads) | 2, 3, 4, 5 | 2, 3, 4, 5 | 2, 3, 4, 5 | 2, 3, 4, 5 | 2 & 4 | 2 & 4 |
| General | | | | | | |
| Safety | Security to open circuit (TTL level) | Security to open circuit (TTL level) | Security to open circuit (TTL level) | Security to open circuit (TTL level) Temperature probe included | Security to open circuit (TTL level) Temperature probe included | Security to open circuit (TTL level) Temperature probe included |

Chassis Specifications

| Essential | SP-50e | SP-150e | VSP | VSP-3e | VMP3e |
|--------------------|--------------------|--------------------|-------------------|--------------------|--------------------|
| Channels available | 1 | 2 | 5 | 8 | 16 |
| Interfaces | Ethernet, USB 2.0 | Ethernet, USB 2.0 | Ethernet, USB 2.0 | Ethernet, USB 2.0 | Ethernet, USB 2.0 |
| Dimension HxWxD | 209 x 136 x 372 mm | 209 x 136 x 372 mm | 95 x 435 x 335 mm | 405 x 225 x 320 mm | 262 x 495 x 465 mm |
| Weight | 3.9 kg | 3.9 kg | 8 kg | 12 kg | 15 kg |
| Power Requirement | 110 W | 110 W | 300 W | 1000 W | 850 W |

A potentiostat for every possible application.

ENERGY STORAGE & CONVERSION

Batteries
Fuel cells & electrolyzers
Supercapacitors
Photovoltaics
Redox Flow Batteries

RESEARCH ELECTROCHEMISTRY

Analytical Electrochemistry
Sensors
Corrosion
MATERIAL SCIENCE



With the largest, most comprehensive range of potentiostats of any manufacturer, you can be sure to find a BioLogic instrument that suits your application.

Here to help.

Online/offline - wherever you are...

BioLogic prides itself in the quality of its potentiostats. We build robust, reliable instruments designed to withstand the rigors of time and the laboratory. But if you do ever encounter a problem with your instrument, you can rest assured that our global support network will be close at hand to help find you a solution quickly and effectively.

And if you just need more information, or perhaps just a little inspiration to help you with your project, you can browse our ever-growing support database of over 500 Learning Centre articles, application/technical notes and support videos at www.biologic.net.

Need high-level EIS measurements? Check out our Premium Range at www.biologic.net

Innovation.

Innovation is engrained in our commercial DNA. The first multi-channel computer-controlled potentiostat (MacPile, 1991), Ethernet connectivity and Embedded EIS are just some of the BioLogic innovations helping scientists around the globe. Our high-quality, high-performance instruments have played a pivotal role in leading research projects since 1983.

www.biologic.net/about us

Shaping the future.
Together.